



# USER MANUAL

PRONTO-50 and PRONTO-250 series | Pocket-Size Power Probes

## **WARRANTY**

### **First Year Warranty**

The Gentec-EO thermal power detectors carry a one-year warranty (from date of shipment) against material and/or workmanship defects when used under normal operating conditions. The warranty does not cover recalibration, or damages related to misuse.

Gentec-EO will repair or replace at our option any PRONTO which proves to be defective during the warranty period; except in the case of product misuse.

Any unauthorized alteration or repair of the product is also not covered by the warranty.

The manufacturer is not liable for consequential damages of any kind.

In the case of a malfunction, contact your local Gentec-EO distributor or the nearest Gentec-EO office to obtain a return authorization number. Return the material to the appropriate address below.

### **Contacting Gentec Electro-Optics Inc.**

To help us answer your calls more efficiently please have the model number of the detector you are using ready before calling Customer Support.

Gentec Electro-Optics, Inc.  
445, St-Jean-Baptiste, Suite 160  
Québec, QC  
Canada, G2E 5N7

Tel: (418) 651-8003  
Fax: (418) 651-1174  
E-mail: [service@gentec-eo.com](mailto:service@gentec-eo.com)  
Website: [gentec-eo.com](http://gentec-eo.com)

## **SAFETY INFORMATION**

Do not use the PRONTO if the device or the detector looks damaged, or if you suspect that the PRONTO is not operating properly.

The user must wait for a while before handling these detectors after power is applied. Surfaces of the detector get very hot and there is a risk of injury if they are not allowed to cool down.

**NOTE:** This equipment has been tested and found to comply with the limits for a Class A digital device, pursuant to part 15 of the FCC Rules. These limits are designed to provide reasonable protection against harmful interference when the equipment is operated in a commercial environment. This equipment generates, uses, and can radiate radio frequency energy and, if not installed and used in accordance with the instruction manual, may cause harmful interference to radio communications. Operation of this equipment in a residential area is likely to cause harmful interference in which case the user will be required to correct the interference at his own expense.

**Caution:** Changes or modifications not expressly approved in writing by Gentec-EO Inc. may void the user's authority to operate this equipment.

## TABLE OF CONTENTS

<b>WARRANTY .....</b>	<b>2</b>
<b>SAFETY INFORMATION .....</b>	<b>3</b>
<b>TABLE OF CONTENTS .....</b>	<b>4</b>
<b>1. PRONTO-50-W5 and PRONTO-250/250-PLUS Portable Laser Probes .....</b>	<b>6</b>
1.1. Introduction.....	6
1.2. Warnings and Disclaimer .....	6
1.3. Specifications .....	7
1.4. Minimum Beam Diameter (Damage Threshold) .....	10
1.5. Mechanical Description .....	11
<b>2. Operating Instructions .....</b>	<b>12</b>
2.1. User Interface.....	12
2.2. Turning the Device ON and OFF .....	12
2.3. Changing the SETTINGS .....	12
2.3.1. Opening and Closing the SETTINGS Menu .....	12
2.3.2. Settings Icons.....	13
2.3.3. Changing the ORIENTATION of the Screen .....	13
2.3.4. Changing the BRIGHTNESS of the Screen.....	13
2.3.5. Adjusting the CORRECTION FACTOR .....	13
2.3.6. Getting Information ABOUT the Device .....	14
2.4. Changing the WAVELENGTH.....	14
2.5. Making a Measurement (SSP mode only) .....	15
2.6. Acquiring, Transferring and Deleting Data .....	16
2.6.1. Acquiring Data .....	16
2.6.2. Transferring Data .....	16
2.6.3. Deleting Data .....	16
<b>3. Operating Instructions (PRONTO-250-PLUS Only).....</b>	<b>17</b>
3.1.1. ZEROING (CWP mode only) .....	17
3.1.2. Setting the SCALE (CWP & SSE modes only) .....	17
3.1.3. Setting the TRIGGER LEVEL (SSE mode only).....	17
3.2. Making a Measurement.....	18
3.2.1. SSP Mode.....	18
3.2.2. CWP mode.....	18
3.2.3. SSE mode .....	19
<b>4. Safety Instructions.....</b>	<b>20</b>
4.1. General.....	20
4.2. Damage to the Optical Absorber Material .....	20
4.3. Not Exceeding the Maximum Sensor Temperature .....	20
4.4. How to Properly Handle the Device .....	21
<b>5. USB Communication .....</b>	<b>22</b>
5.1. Description .....	22
5.2. Setting up communication to the PRONTO .....	22
5.2.1. Verify the COM Port.....	22
5.2.2. Connect the PRONTO .....	22
5.2.3. To Echo Commands .....	23
5.2.4. Test the Connection .....	23
5.2.5. To Add a Character Delay .....	23
5.2.6. HyperTerminal Settings Shortcut.....	23
5.3. Serial Command Format .....	23
5.3.1. Serial Protocol Rules .....	23

5.3.2.	Text Mode Rules .....	23
5.4.	List of Serial Commands for the PRONTO (Summary) .....	24
5.5.	Detailed Description of the Serial Commands for PRONTO (Complete).....	25
5.5.1.	Display .....	25
5.5.2.	Data Acquisition .....	29
5.5.3.	Setup.....	31
5.5.4.	Control.....	32
5.5.5.	Instrument and Detector Information .....	37
5.5.6.	Error Messages.....	41
<b>6.</b>	<b>USB Driver Installation .....</b>	<b>42</b>
<b>7.</b>	<b>Maintenance .....</b>	<b>42</b>
7.1.	Free Firmware Upgrade .....	42
<b>8.</b>	<b>Declaration of Conformity .....</b>	<b>43</b>
<b>9.</b>	<b>UKCA Declaration of Conformity.....</b>	<b>44</b>
	<b>Appendix A: WEEE Directive .....</b>	<b>45</b>
	<b>Appendix B: Installing the PRONTODataTransfer Software .....</b>	<b>46</b>
	<b>Appendix C: Updating the PRONTO Firmware.....</b>	<b>48</b>

## 1. PRONTO-50-W5 AND PRONTO-250/250-PLUS PORTABLE LASER PROBES

### 1.1. INTRODUCTION

The PRONTO-50-W5 and PRONTO-250 Portable Laser Probes from Gentec-EO are pocket-size power and energy meters with built-in display. The PRONTO-50-W5 and PRONTO-250 give accurate measurement of laser power up to 50 W and 250 W in only 5 seconds. In addition to the regular Single Shot Power (SSP) mode, The PRONTO-250-PLUS features 2 additional measurement modes: in Continuous Power (CWP) mode, the device measures power continuously up to 8 W while in Single Shot Energy Mode (SSE), it measures the energy of individual pulses. The high power surface absorber sensors are designed to be used at high average power densities. PRONTO detectors can be supplied with an optional stand and/or power supply.

If the desired wavelength is outside the calibrated spectral range, you can use the “*Correction Factor*” function to adjust the displayed measures.

Call your nearest Gentec-EO distributor to repair and/or to recalibrate the PRONTO. To find the nearest Gentec-EO office or distributor in your country, refer to p. II, **Contacting Gentec Electro-Optics Inc.**

### 1.2. WARNINGS AND DISCLAIMER

Laser power detection is completely automatic. There is no need for an external timer.



#### Warnings

WHEN **HOT** APPEARS ON THE SCREEN, REMOVE THE DEVICE FROM THE BEAM IMMEDIATELY TO AVOID DAMAGING THE SENSOR.

Be careful not to exceed the maximum levels and densities stated in the specifications.

The handle on the PRONTO is for carrying purposes only. The user must not hold the PRONTO during measurements. Extreme caution must be taken during laser power measurements.

In no event shall Gentec-EO or any of its affiliates be liable for any indirect, special, incidental or consequential injury to persons or damage to property caused by the use of any of our products, by purchasing from Gentec-EO or any of its affiliates, you hereby indicate that you understand and agree to the following:



#### Disclaimer

I am fully responsible for the safe application and use of this detector and agreed to such by completing the sales process.

I will not use a laser device without wearing approved laser safety goggles designed for such purpose.

I am aware and responsible of safely dealing with any back reflections.  
I will not use the detector in violation of any local, state or federal law, and  
I understand that it is my responsibility to know and abide by those laws relating to the ownership and use of the detector in my jurisdiction.

### 1.3. SPECIFICATIONS

These products all have a calibration at 1064 nm (YAG) and they also have a traceable wavelength correction to complete the calibrated spectral range. The PRONTO-250/250-PLUS have an additional calibration at 10.6  $\mu\text{m}$  ( $\text{CO}_2$ ). The PRONTO-250-PLUS units also have an additional calibration for SSE mode.

The following specifications are based on a one-year calibration cycle, an operating temperature of 15 to 28°C (59 to 82°F) and a relative humidity not exceeding 80%. Monitors must be stored in an environment between 10°C to 60°C (50 to 140°F) and a relative humidity not exceeding 90%.

	PRONTO-250	PRONTO-250-PLUS	PRONTO-50-W5
	<b>Sensor &amp; Calibration</b>		
<b>Aperture</b>	19 mm Ø		
<b>Spectral Range</b>	0.19 - 20 $\mu\text{m}$		0.19 - 10 $\mu\text{m}$
<b>Calibrated Spectral Range<sup>1</sup></b>	0.248 - 2.5 $\mu\text{m}$ and 10.6 $\mu\text{m}$		0.248 - 2.5 $\mu\text{m}$
<b>Programmed Spectral Range<sup>2</sup></b>	0.193 - 10.6 $\mu\text{m}$		0.193 - 2.5 $\mu\text{m}$
<b>Power Range (in SSP mode)</b>	0.5 - 250 W		0.5 - 50 W
<b>Noise Level</b>	10 mW		4 mW
<b>Response Time (in SSP mode)</b>	5 s		
<b>Calibration Uncertainty</b>	$\pm 3 \%$		
	<b>Damage Thresholds &amp; Laser Limits</b>		
<b>Maximum Average Power Density</b> 1064 nm, 10 W, CW 10.6 $\mu\text{m}$ , 10 W, CW	45 kW/cm <sup>2</sup> 14 kW/cm <sup>2</sup>		100 kW/cm <sup>2</sup>
<b>Maximum Number of Readings and Exposure Time Before Probe Must Be Cooled</b> (for 25°C starting temp. see section 4.3 for more information)	<b>Power</b>	<b>Max # Readings / Max Exposure Time</b>	
	8 W	$\infty$	$\infty$
	25 W	10	60 sec
	50 W	5	30 sec
	125 W	2	12 sec
	250 W	1	6 sec
<b>Maximum Device Temperature at Maximum Power</b>	65°C		
<b>Pulsed Laser Damage Thresholds : Maximum Energy Density</b> 1064 nm, 360 $\mu\text{s}$ , 5 Hz 1064 nm, 7 ns, 10 Hz 532 nm, 7 ns, 10 Hz 266 nm, 7 ns, 10 Hz	9 J/cm <sup>2</sup> 1.0 J/cm <sup>2</sup> 0.6 J/cm <sup>2</sup> 0.3 J/cm <sup>2</sup>		100 J/cm <sup>2</sup> 1.1 J/cm <sup>2</sup> 1.1 J/cm <sup>2</sup> 0.7 J/cm <sup>2</sup>
<b>Pulsed Laser Damage Thresholds : Peak Power Density</b>  1064 nm, 360 $\mu\text{s}$ , 5 Hz 1064 nm, 7 ns, 10 Hz 532 nm, 7 ns, 10 Hz 266 nm, 7 ns, 10 Hz	25 kW/cm <sup>2</sup> 143 MW/cm <sup>2</sup> 86 MW/cm <sup>2</sup> 43 MW/cm <sup>2</sup>		667 kW/cm <sup>2</sup> 157 MW/cm <sup>2</sup> 157 MW/cm <sup>2</sup> 27 MW/cm <sup>2</sup>
<b>Minimum Frequency for Pulsed Lasers</b>	15 Hz		

<sup>1</sup> Pronto-50-W5 standard units are calibrated at one wavelengths (1064nm) and have a NIST-traceable wavelength correction factor to cover the complete calibrated spectral range.  
Pronto-250 standard units are calibrated at two wavelengths (1064nm and 10.6 $\mu\text{m}$ ) and have a NIST-traceable wavelength correction factor to cover the complete calibrated spectral range.  
In addition to the standard Pronto-250 calibrations, the Pronto-250-PLUS standard units are also calibrated for SSE measurements.

<sup>2</sup> Available wavelengths in the user interface.

	Physical Characteristics
Dimensions	59W x 181L x 17D mm
Weight	210 g
Mounting Holes (for post)	2 x 8-32
Operating Conditions	Ambient temperature: 15-28°C Maximum relative humidity: 80%
Storage Conditions	Ambient temperature: 10-60°C Maximum relative humidity: 90%
Battery Type	USB rechargeable Li-ion
Battery Life	17 hours or 4200 measurements (with brightness set at 25%) Charge time: 7.5 hours when totally empty
Maximum Battery Cycle Count (Number of charge cycles before the battery loses its efficiency)	Approximatively 500 full charges (0-100%)
Lithium Battery Transport Compliance	UN: Recommendations on the Transport of Dangerous Goods, Manual of Tests and Criteria, ST/SG/AC.10/11/Rev.6/Amend.1 Section 38.3 <sup>1</sup>
	User Interface
Interface	Touchscreen
Display	Color LCD 28 X 35 mm (128 x 160 pixels)
Data Acquisition and Transfer	Maximum of 50 000 measurements Acquisition frequency of 6.8Hz in CWP mode
Screen Personalization	4 screen orientations and 4 brightness levels
Saved Settings	Screen Orientation, Screen Brightness, Wavelength and Correction Factor
Display Numerical Resolution	4 digits

	PRONTO-250-PLUS
	CWP Mode
Power Range in CWP mode	0.01 - 8 W
Noise Level in CWP mode	10 mW (Excluding possible offset and used with an opto-mechanical stand)
Typical Rise Time in CWP mode (0-95%)	1.5 s
Maximum Device Temperature at Maximum Power (in CWP mode)	55°C
Calibration Uncertainty in CWP mode	± 2.5 %
	SSE Mode
Energy Range in SSE mode (>1 ns)	0.06 - 25J up to 150J for pulse >1ms
Noise Level in SSE mode	60 mJ
Trigger Level	0.1 - 99.9 % (Default level is 2%)
Minimum Repetition Period	4 s
Maximum Pulse Width	88 ms
Typical Rise Time in SSE mode (0-100%)	0.26 s
Calibration Uncertainty in SSE mode	± 5 %

Specifications are subject to change without notice


<sup>1</sup> UN38.3 certificate available upon request.






#### 1.4. MINIMUM BEAM DIAMETER (DAMAGE THRESHOLD)

The following graphic shows the diameter corresponding to the damage threshold for a Gaussian beam profile. The “minimum  $1/e^2$  beam diameter” is calculated to obtain a peak intensity 50% lower than the damage threshold and should be considered as the “safe” minimum diameter.

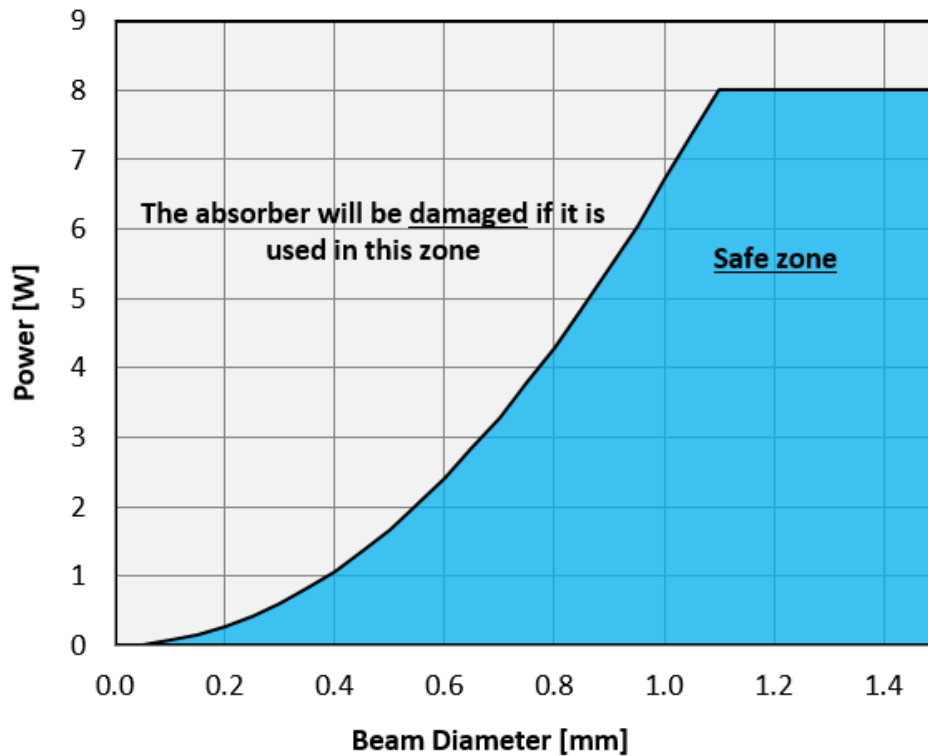


**Figure 1** Minimum Beam Size for PRONTO-250-PLUS absorber at 1064 nm in CWP mode

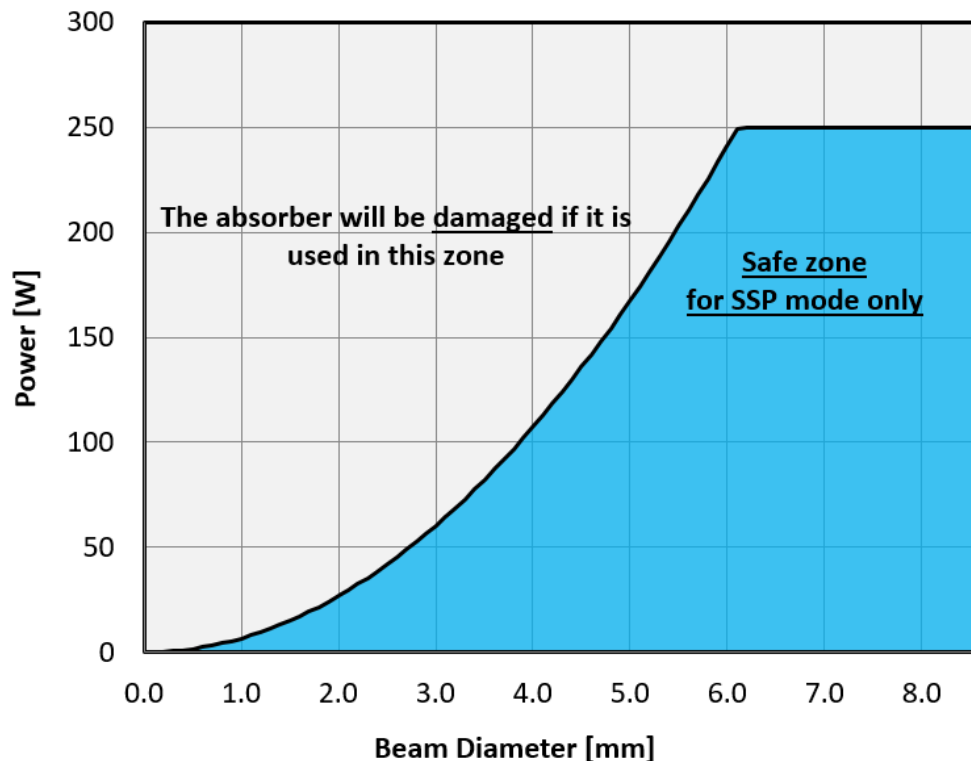

**Figure 2** Minimum Beam Size for PRONTO-250 and PRONTO-250-PLUS absorber at 1064 nm in SSP mode

## 1.5. MECHANICAL DESCRIPTION

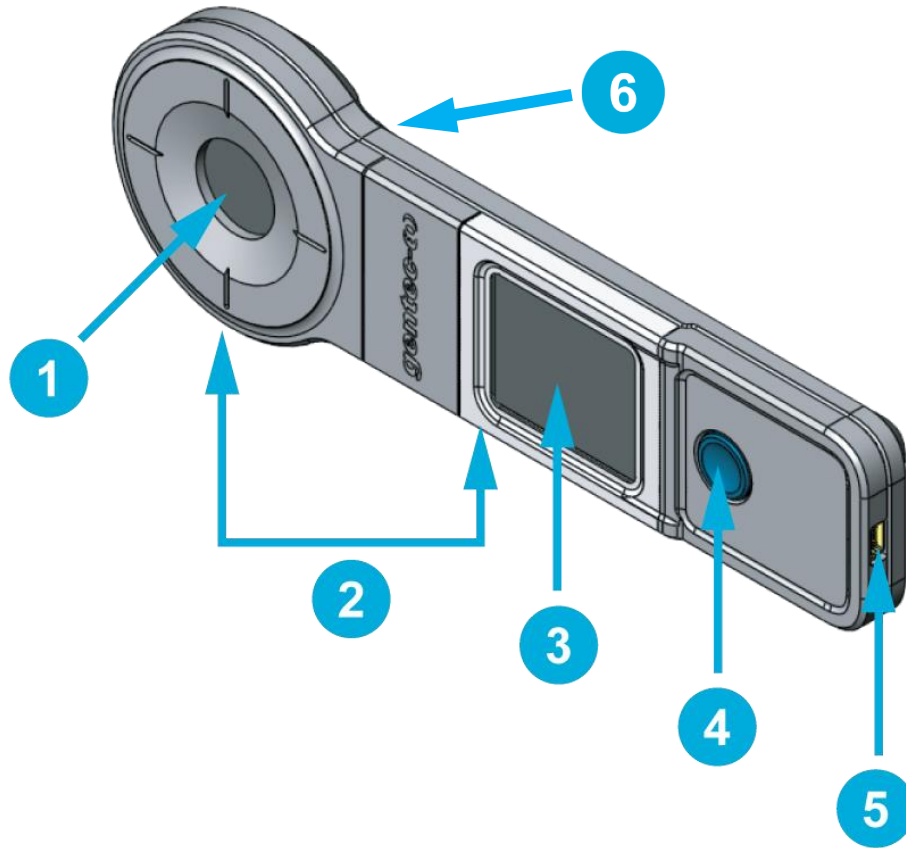

**Figure 3** PRONTO mechanics

**1**

**Absorber**

The laser must be centered on the absorber when making a measurement.

**2**

**Mounting Holes**

There are 2 8-32 mounting holes to fit the device on a post for a safe use during the measurements.

**3**

**Touch Screen Display and Controls**

The touch screen interface controls the device.

**4**

**On/Off/Settings Button**

This button is used to power the device on (press when the device is off) and off (press and hold for 3 seconds when the device is on). It is also used to access the Settings menu (press when the device is on).

**5**

**USB Port**

The Mini-B USB2.0 port is used to transfer data from the device to a PC and to charge the battery.

**6**

**Overheating Indicator**

The central part of the yellow sticker on the rear face turns to black when the temperature of the device exceeds 65 °C (150 °F). This indicates that the detector was probably damaged due to overheating. In this case, contact your local Gentec-EO representative.

## 2. OPERATING INSTRUCTIONS

### 2.1. USER INTERFACE

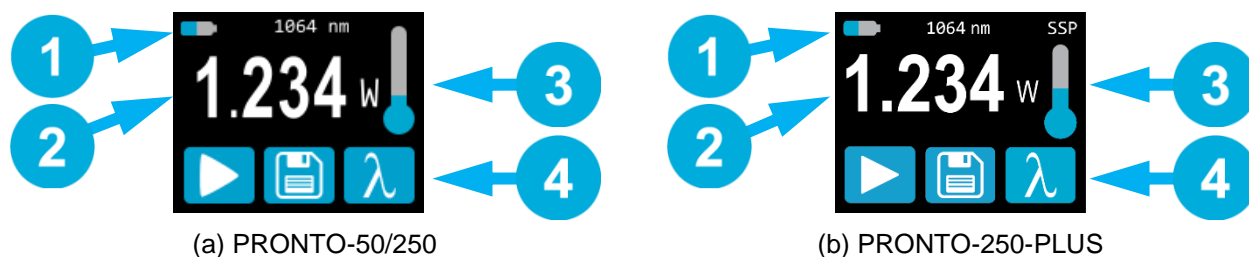


Figure 4 User Interface

1

#### Device State

The top portion of the display shows the battery level and selected wavelength. In the PRONTO-250-PLUS, the top portion also indicates the measurement mode. Its default mode is SSP, which works like the PRONTO-50/250.

2

#### Measured Value

The central part of the display shows the last measured value. In the PRONTO-250-PLUS, this section also serves as a button to access the scales menu in CWP and SSE modes.

3

#### Temperature

The thermometer icon shows the head temperature.



*If the thermometer is red and the screen displays **HOT**, block the laser IMMEDIATELY and let the instrument cool down.*

4

#### Controls

The first button at the bottom of the display is specific to the measurement mode. The second button (Save) is for data acquisition, and the third button (Wavelength) opens the Wavelength menu.

### 2.2. TURNING THE DEVICE ON AND OFF

**Turn ON:** Press the **On/Off/Settings** button



**Turn OFF:** Press and Hold the **On/Off/Settings** button for 3 seconds




**Tip**

The device will automatically turn off after 5 minutes of inactivity, except if data is being acquired. When the device is plugged into a PC via the USB cable, it will power on and stay on until it is manually turned off or unplugged.

### 2.3. CHANGING THE SETTINGS

#### 2.3.1. Opening and Closing the SETTINGS Menu

When the device is on, short-press the **On/Off/Settings** button  to access the menu. Pressing the button again will close it and save whatever settings have been changed.

### 2.3.2.Settings Icons

The following icons are presented in the Settings menu. The measurement mode options are only available with the PRONTO-250-PLUS, since PRONTO-250 and PRONTO-50-W5 always works in Single-Shot Power mode.

ICON	NAME	DESCRIPTION
	Screen Orientation	Indicates the <b>Screen Orientation</b> (4 options) Press → Toggles to the next orientation
	Brightness	Indicates the screen's <b>Brightness</b> level (4 options) Press → Toggles to the next brightness level
	Correction Factor	Press → Opens the <b>Correction Factor</b> (User Calibration) menu
	About	Press → Opens the <b>About</b> menu
FOR PRONTO-250-PLUS ONLY		
	More	Press → Switches to the secondary menu
	Single Shot Power	Press → Goes to <b>Single Shot Power</b> Mode
	Continuous Power	Press → Goes to <b>Continuous Power</b> Mode
	Single Shot Energy	Press → Goes to <b>Single Shot Energy</b> Mode

### 2.3.3.Changing the ORIENTATION of the Screen

There are 4 possible screen orientations: Up, Down, Left and Right. The button displays the current orientation, for example Up: . To change the orientation, simply press the button and the next choice will appear. The changes will be saved once you exit the Settings menu.


### 2.3.4.Changing the BRIGHTNESS of the Screen

There are 4 possible brightness levels: 100%, 75%, 50% and 25%. The button displays the current brightness, for example 75%: . To change the brightness, simply press the button and the next choice will appear. The changes will be saved once you exit the Settings menu.

### 2.3.5.Adjusting the CORRECTION FACTOR


It is possible to adjust the calibration of the device. You can do so by applying a **Correction Factor** to the measurements. To access the **Correction Factor** menu, simply press the CAL button . The number displayed is the multiplication factor that will be applied to the measurements. A factor of 1.000 keeps the calibration unchanged. You can enter any number between 0.500 and 2.500.


The 4 digits of the correction factor are changed individually. For example, if the **Correction Factor** is 1.000 and you want to change it to 1.200, just press the first 0 (make sure it turns blue, like this ) and then increment it twice with the Up button.

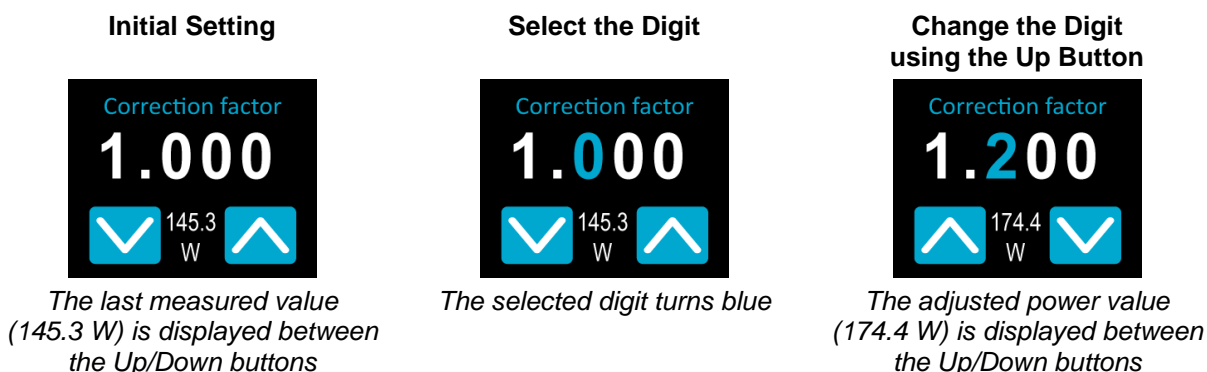

Once the value is entered, press the On/Off/Settings button to exit the menu and automatically save your new settings.

### 2.3.6. Getting Information ABOUT the Device


Relevant information about the device are stored in the **About** menu . This information can be useful when returning the device for repairs or recalibration. These include, the **Model Name**, the **Serial Number**, the **Firmware Version** and the **Last Calibration Date**.

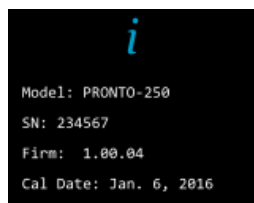

**Figure 5** The About menu shows information about the device

Press the On/Off/Settings button to exit the menu and go back to the measurement screen.



## 2.4. CHANGING THE WAVELENGTH

Before making a measurement, it is important that you adjust the calibration of the device to the wavelength of the source to be measured. To do so, press the **Wavelength** button  to display a list of presets.



**Figure 6** The 6 wavelength buttons can be customized by the user

There are 6 presets in the wavelength menu. The values can be set to any wavelength defined by the programmed spectral range. To change a value, press it until the screen changes to the **Set Wavelength** menu. Each digit is changed individually by selecting it and pressing the Up/Down arrows. For example, if you want to change 10.6 μm to 532 nm, you need to do the following:

1. Select the units. Make sure they turn blue  and use the Up or Down button to change them to nm
2. Select the 1st number. Make sure it turns blue  and use the Down button to reduce it to 0

3. Select the 2nd number. Make sure it turns blue **0060nm** and use the Up button to increase it to 5
4. Select the 3rd number. Make sure it turns blue **0560nm** and use the Down button to reduce it to 3
5. Select the 4th number. Make sure it turns blue **0530nm** and use the Down button to reduce it to 2
6. You're done! **0532nm**

Once the desired value is entered, press the On/Off/Settings button to exit the menu and automatically save your new settings.

## 2.5. MAKING A MEASUREMENT (SSP MODE ONLY)


Once all the settings are adjusted, you are ready to make a measurement. Just follow the step-by-step instructions below.




### Tip

We recommend placing the device on an optical stand or on a horizontal surface when making a measurement.


Be careful to conform to all the recommended specifications for beam size, placement and laser power.

1. Turn the device ON.
2. Press the PLAY  button, a sequence of dots will appear. This indicates that the device is waiting for a laser beam to trigger the measurement.




**NOTE:** At any time during the measurement sequence, you can press the STOP  button. The device will stop the measurement process and go back to its initial state (displaying the last measured value).

3. Place the device in the laser beam path, with the laser beam centered on the absorber.
4. As soon as the device detects a laser beam, it will automatically start the 5 seconds countdown for the measurement. Leave the device in the beam path for the entire countdown period, which will be indicated on the screen.

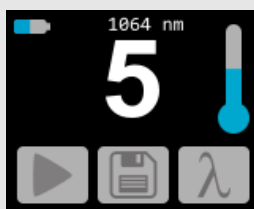

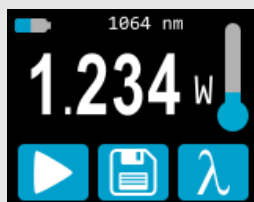


5. Once the countdown is finished, the measurement will appear. You can remove the device from the beam path.



6. The reading will stay on the display until the next measurement, even if you turn the device off and on again.
7. To make a new measurement, go back to Step 2.


## 2.6. ACQUIRING, TRANSFERRING AND DELETING DATA

### 2.6.1.Acquiring Data

You can store the measurements done by the device simply by pressing the **Save** button . When pressed, the button turns white  to indicate that data is being stored in the internal memory of the device. Once activated, the data acquisition will stay active until stopped. To stop the data acquisition, press the Save button again and it will revert to its original state .

### 2.6.2.Transferring Data

To retrieve the data, you must connect the PRONTO to your computer with a USB cable and use the

**PRONTODataTransfer** software . You can download our latest version of the software in the *Downloads* section of our website (<https://gentec-eo.com/downloads>). The data will be uploaded on your computer in a text format, which you can save to a known location on your computer and then open in your preferred analysis software.




#### Warning

Once data has been transferred to a computer, it is deleted from the internal memory of the device.

Complete installation and data transfer instructions can be found in **Appendix B: Installing the PRONTODataTransfer Software**.


### 2.6.3.Deleting Data


To delete data from the internal memory, you must hold the **Save** button  approximately 3 seconds. A text will appear and ask **Erase all data**. Hit the **YES** button than click **OK**.




### 3. OPERATING INSTRUCTIONS (PRONTO-250-PLUS ONLY)

#### 3.1.1.ZEROING (CWP mode only)

1. Block off any laser radiation to the detector.
2. To set the zero, wait until the reading has stabilized. The power read by the device when no laser beam is incident on the sensor may not be exactly zero if the device is not thermally stabilized. Wait until the reading without laser power is stable for several minutes. Half an hour warm-up is recommended for measuring low powers precisely.
3. Press the ZERO button . While the device is busy, the control buttons are inactive.

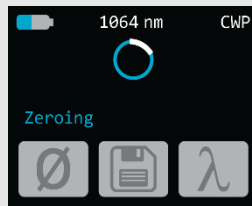

4. When the process is complete, the device returns to the measurement screen and you are now ready to make an accurate measurement.

When the offset is applied to the measurement, the **Zero** button is white . The offset stays active until the **Zero** button is pressed again or until the PRONTO is turned off.

#### 3.1.2.Setting the SCALE (CWP & SSE modes only)

To access the **Range** menu, you must press on the measured value (the central part of the touchscreen) for approximately 3 seconds. The Up/Down arrows scroll through the available ranges and AUTO, which is the default, automatic range setting.




**Figure 7** In the **Range** menu, use the Up and Down buttons to set the manual range or to go back to the default “Auto” range setting.




Once the desired value is entered, press the On/Off/Settings button to exit the menu and automatically save your new settings.

#### 3.1.3.Setting the TRIGGER LEVEL (SSE mode only)

In SSE mode, it is possible to control the trigger level. To access the **Trigger Level** menu, simply press the

**Trigger Level** button  at the bottom of the measurement screen. Accepted values are between 0.1% and 99.9%. The default value is 2.0%.

The 3 digits of the trigger level are changed individually. For example, if the **Trigger Level** is 2.0% and you want to change it to 5.0%, just press the 2 (make sure it turns blue) and then increment it three times with the Up button.

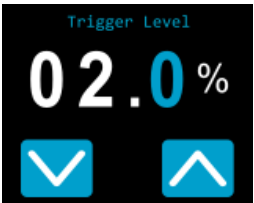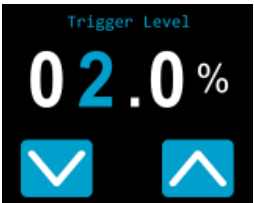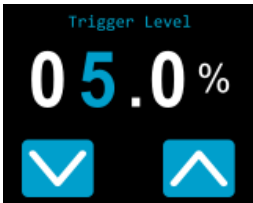Initial Setting	Select the Digit	Change the Digit using the Arrows
		
<i>The default value is 2.0%</i>	<i>The selected digit turns blue</i>	<i>The adjusted level is saved automatically when exiting this screen</i>

Once the desired value is entered, press the On/Off/Settings button to exit the menu and automatically save your new settings.

### 3.2. MAKING A MEASUREMENT

Once all the settings are adjusted and the desired measurement mode is selected, you are ready to make a measurement. Just follow the step-by-step instructions below.



#### Tip

We recommend placing the device on an optical stand or on a horizontal surface when making a measurement.


Be careful to conform to all the recommended specifications for beam size, placement and laser power.

#### 3.2.1.SSP Mode

Instructions for this mode are detailed in Section 2.5.



#### 3.2.2.CWP mode

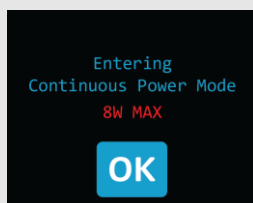1. Turn the device ON.
2. The PRONTO displays a warning (8W maximum), press OK.



3. For maximum accuracy, adjust the reading to zero before making measurements (see Section 3.1.1). For low power measurements, the sensor must be protected from ambient light and temperature fluctuations.
4. Place the device in the laser beam path, with the laser beam centered on the absorber.
5. The device will automatically display and update the power value.

### 3.2.3.SSE mode

1. Turn the device ON.
2. Place the device in the laser beam path, with the laser beam centered on the absorber.
3. As soon as the device detects a laser pulse, it will automatically measure its energy.
4. The TRIGGER LEVEL button  turns gray  every time a pulse is detected.
5. For accurate results, do not touch the screen while taking measurements and conform to the specifications for laser pulse width and repetition rate.

## 4. SAFETY INSTRUCTIONS

### 4.1. GENERAL

To ensure a long lifetime of accurate measurements, the PRONTO power probe should be maintained within the following ambient conditions:

Storage environment temperature: 10 to 60°C, RH < 90%  
Operating environment temperature: 15 to 28°C, RH < 80%.

It is possible to store and operate your PRONTO power probe beyond this range. For any specific requirements, please contact your local Gentec-EO representative.

For the most accurate measurements, center the beam on the sensor. Ideally, the beam diameter should be the same as the one used during calibration. The beam diameter at calibration corresponds to >98% of the encircled power centered on 50% of the sensor's surface area. This complies with the International Electrotechnical Commission standard #1040: "Power and Energy Measuring Detector [...]". Refer to the calibration certificate for the exact beam diameter used during calibration.

### 4.2. DAMAGE TO THE OPTICAL ABSORBER MATERIAL

Damage to the optical absorber material is usually caused by exceeding the manufacturer's specified maximum incident **Average Power Density**. Refer to the specifications table.

The PRONTO-250-PLUS series can measure up to 250W. The beam diameter should always be as large as possible to avoid damaging the absorber. **We recommend using a beam size diameter that covers 70% to 80% of the sensor's aperture.** For the PRONTO, this corresponds to a diameter of 16-17 mm.

In any case, the beam's incident area should not be less than 10% of the detector's area. Please contact Gentec-EO to make measurements with smaller beams.

The damage threshold decreases with the laser power. Refer to the graphic in Figure 2 for the diameter corresponding to the damage threshold for a Gaussian beam profile. The "minimum 1/e<sup>2</sup> beam diameter" is calculated to obtain a peak intensity 50% lower than the damage threshold and should be considered as the "safe" minimum diameter. If there are "hot spots" in the beam profile, they must be considered in the calculation of the peak intensity.

Damage may also be caused if you use a detector with a contaminated absorber surface. Slight discoloration of the coating may occur, but this does not affect the calibration.

### 4.3. NOT EXCEEDING THE MAXIMUM SENSOR TEMPERATURE

The PRONTO has a built-in thermometer to monitor the sensor's temperature. In order to avoid damaging the sensor, its temperature must not exceed the maximum allowable absorber temperature as in the specification. For temperatures of 25°C or lower, the thermometer's bulb will be filled as shown in the picture at the right of this text. When making a measurement, the sensor's temperature will gradually rise and the thermometer icon will fill up too.



If the temperature of the sensor gets close to the acceptable limit, the screen will display the mention HOT in bold red letters, the thermometer icon will be red and all functions will be disabled.



*If this screen appears, it is essential that you remove the device from the laser beam path and allow it sufficient time to cool before attempting another measurement.*



In addition to the built-in thermometer, the PRONTO has a yellow temperature sticker on its rear face. The central part of the sticker turns to black when the temperature of the device exceeds 65 °C (150 °F). This indicates that the detector was probably damaged due to overheating. In this case, contact your local Gentec-EO representative.

#### 4.4. HOW TO PROPERLY HANDLE THE DEVICE

Laser radiation can heat the device very quickly. The entire body of the device can get hot, especially at higher powers. In addition to the standard safety precautions that should be taken when working with laser radiation, we recommend placing the device on an optical stand when making a measurement. For this use, the device is equipped with 2 mounting holes with an 8-32 thread (see Figure 3).



**Figure 8** Recommended setup when making a measurement: install the device on an optical stand before placing it in the laser beam path.

Gentec-EO offers a compatible stand for the PRONTO, you can contact your local Gentec-EO representative for the exact model.

It is also possible to place the PRONTO on a horizontal surface and illuminate the sensor from above. The back of the device is flat so it will stay stable when placed on a horizontal surface.



**Figure 9** Recommended setup when making a measurement: install the device flat on a horizontal surface and illuminate from above.

## 5. USB COMMUNICATION

### 5.1. DESCRIPTION

The PRONTO has one communication mode and it is in ASCII. The format will require text input commands which follow rules stated in section 5.3. The section 5.5 describes all the commands.

The USB class used by the PRONTO is a CDC, or Communications Device Class. This means it shows up in the host PC as a COM port, but it is not a COM port, rather a true full speed USB port. You can talk to it as if it were an RS232 port, but much faster. Follow the Windows Prompts to install the USB drivers. The USB drivers are fully tested and digitally signed by Microsoft.

Open the appropriate port in your software with standard COM port tools. None of the port settings matter since they are not used, so leave them at whatever default they are in. It's a real USB connection.

Use the standard COM port write and read protocols to control the PRONTO.

### 5.2. SETTING UP COMMUNICATION TO THE PRONTO

#### 5.2.1. Verify the COM Port

To verify the USB installation and find the COM port number, click:

**Start → Settings → Control Panel → System → Device Manager**

Scroll down to **Ports (COM & LPT)** and double click that line. One of the options should be

**Power/Energy Monitor (COM#)**

Note the COM port number, you will need it for the next step.

#### 5.2.2. Connect the PRONTO

You may use any serial communications software that you are familiar with. Our instructions are for HyperTerminal because it is widely available on PCs with Windows™.

Select:

**Start → Programs → Accessories → Communications → HyperTerminal**

To save communication settings, enter a name for the connection. In the drop-down menu for **“Connect using”** select the COM port that the USB driver was installed on (Section 5.2.1). Select **OK**.

Input the following settings into the communications parameter window that appears next.

PRONTO COM Port Settings	
Bits per second	Any settings will work
Data bits	Any settings will work
Parity	Any settings will work
Stop bits	Any settings will work
Flow control	Any settings will work

Click **OK** to begin entering serial commands in the HyperTerminal window.

### 5.2.3. To Echo Commands

The commands you type will not appear in the HyperTerminal window unless you set up HyperTerminal to do so. Only the response from the PRONTO will be displayed. If you prefer to see the commands you are typing on the HyperTerminal window click the **File** menu and execute the following sequence:

**File → Properties → Settings → ASCII setup → “Echo typed characters locally” → OK**

### 5.2.4. Test the Connection

In the HyperTerminal window, type **\*VER**. If the response is the version of your PRONTO, you are successfully connected and ready for serial command action.

### 5.2.5. To Add a Character Delay

In some cases, the PRONTO has connection problems that can be solved by adding a character delay. To adjust this delay, click the File menu and execute the following sequence:

**File → Properties → Settings → ASCII setup → “Character delay”: add a few milliseconds of delay → OK**

### 5.2.6. HyperTerminal Settings Shortcut

When you end the session, HyperTerminal asks if you want to save your settings. To avoid inputting the communication parameters again in the future, save by clicking **Yes**. The next time you execute the string of commands, the name of your session will appear after HyperTerminal. Clicking on the session name will open the connection using the saved settings. To avoid re-entering the string of commands, put a shortcut to this file on your desktop: search for the file name, select the file. Right click and select **Shortcut** in the drop down menu.

## 5.3. SERIAL COMMAND FORMAT

### 5.3.1. Serial Protocol Rules

Commands are sent as text strings. The response will either be data or an empty string.

### 5.3.2. Text Mode Rules

All text commands must begin with a trig character (\*). You do not need to end with a line-feed and/or a carriage-return. Parameters must NOT be separated by spaces. Characters do not have to be capitals, mixed upper and lower cases is ok. Replies to all text mode commands are also in text mode, and end with a carriage-return and a line-feed.

In case of an error, the reply string is one of the following :

**“Command Error. Command not recognized.”**

**or**

**“Command Error. Command must start with '\*'”**

Because all Text Mode replies end with a carriage return <CR> or line-feed <LF> (or both), a text reply contains tabulations when many elements need to be separated in the string. This is useful when exporting data to a spreadsheet.

**5.4. LIST OF SERIAL COMMANDS FOR THE PRONTO (SUMMARY)**

#	Command Name	Command	Description
<b>DISPLAY</b>			
01	Set Scale	SCS	Manually sets the scale
02	Set Scale Up	SSU	Changes scale to the next higher scale
03	Set Scale Down	SSD	Changes scale to the next lower scale
04	Get Current Scale Index	GCR	Returns scale index between 0 and 41
05	Set Autoscale	SAS	Sets the autoscale
06	Get Autoscale	GAS	Returns autoscale status
07	Display Valid Scale	DVS	Displays the valid scales for the connected head
08	Set Trigger Level	STL	Sets the internal trigger level for pulse energy
09	Get Trigger Level	GTL	Returns trigger level value
10	Get Measure Mode Display	GMD	Returns the current measure mode on PRONTO
11	Control LCD	LCD	Turn On/Off the LCD
<b>MEASUREMENT</b>			
<b>Data Acquisition</b>			
12	Query Current Value	CVU	Gets the value currently in ASCII or binary
13	Send Continuous Transmission of Data	CAU	Sends the values in ASCII or binary to the serial port with the data sampling setting
14	Stop the CAU Command	CSU	Stops the *CAU command
15	Query New Value Ready	NVU	Determine if new reading is available or not
16	Get Housing Temperature	TMP	Return the housing temperature
<b>Setup</b>			
17	Set Personal Wavelength Correction in nm	PWC	Specifies the wavelength in nm
18	Set Personal Wavelength Correction in microns	PWM	Specifies the wavelength in microns
19	Get Wavelength	GWL	Returns the wavelength in nm
<b>Control</b>			
20	Set Anticipation	ANT	Turns the anticipation on or off
21	Get Anticipation Status	GAN	Returns the anticipation status
22	Set Zero Offset	SOU	Zeroes the reading for a value without offset
23	Clear Zero Offset	COU	Undoes the zeroing of the reading for a power detector
24	Get Zero Offset	GZO	Returns the zero offset status
25	Set User Multiplier	MUL	Sets the multiplier value
26	Get User Multiplier	GUM	Returns the current multiplier value
27	Set User Offset	OFF	Sets the offset value
28	Get User Offset	GUO	Returns the current offset value
29	Short Touch on Button	PRS	Simulate a short touch on button of display
30	Long Touch on Button	PRL	Simulate a long touch on button of display
<b>INSTRUMENT AND DETECTOR INFORMATION</b>			
31	Query Version	VER	Gets firmware version of the monitor
32	Query Status	STS	Retrieves the detector information and monitor settings
33	Query Extended Status	ST2	Returns the extended status
34	Return Instrument ID	IDN	Returns the device model
35	Return Instrument Firmware	GSV	Returns the firmware version
36	Return Global Information	GFW	Returns the firmware identification number, the device model and firmware version.
37	Query Battery State	QSO	Return the battery level



## 5.5. DETAILED DESCRIPTION OF THE SERIAL COMMANDS FOR PRONTO (COMPLETE)

### 5.5.1. Display

#### 01 - Set Scale

This command is used to force the display of the current data into a specific scale. The lower scale is always zero, the higher scales can be found in the table below. The Autoscale mode applies the best scale for the current values in real time. The parameter must be one of the identifiers in the table below and have 2 digits.

Command	Parameters	Answer	Model Available
SCS	Range index		PRONTO-250-PLUS

Range Identifiers:

Index	Value	Index	Value
00	1 picowatt or picojoule	21	30 milliwatts or millijoules
01	3 picowatts or picojoules	22	100 milliwatts or millijoules
02	10 picowatts or picojoules	23	300 milliwatts or millijoules
03	30 picowatts or picojoules	24	1 watt or joule
04	100 picowatts or picojoules	25	3 watts or joules
05	300 picowatts or picojoules	26	10 watts or joules
06	1 nanowatt or nanojoule	27	30 watts or joules
07	3 nanowatts or nanojoules	28	100 watts or joules
08	10 nanowatts or nanojoules	29	300 watts or joules
09	30 nanowatts or nanojoules	30	1 kilowatt or kilojoule
10	100 nanowatts or nanojoules	31	3 kilowatts or kilojoules
11	300 nanowatts or nanojoules	32	10 kilowatts or kilojoules
12	1 microwatt or microjoule	33	30 kilowatts or kilojoules
13	3 microwatts or microjoules	34	100 kilowatts or kilojoules
14	10 microwatts or microjoules	35	300 kilowatts or kilojoules
15	30 microwatts or microjoules	36	1 megawatt or megajoule
16	100 microwatts or microjoules	37	3 megawatts or megajoules
17	300 microwatts or microjoules	38	10 megawatts or megajoules
18	1 milliwatt or millijoule	39	30 megawatts or megajoules
19	3 milliwatts or millijoules	40	100 megawatts or megajoules
20	10 milliwatts or millijoules	41	300 megawatts or megajoules

**Default:** Autoscale.



#### Example

The following example sets the scale to 30 watts or joules:

Command: *SCS27	Answer:
-----------------	---------

02 - Set Scale Up

This command is used to force the display of the current data into a higher scale.

Command	Parameters	Answer	Model Available
SSU	None		PRONTO-250-PLUS

03 - Set Scale Down

This command is used to force the display of the current data into a lower scale.

Command	Parameters	Answer	Model Available
SSD	None		PRONTO-250-PLUS

04 - Get Current Scale Index

This command returns the scale index between 0 and 41. Please refer to Set Scale command (SCS) details for the complete scale index table.

Command	Parameters	Answer	Model Available
GCR	None	Index from 0 to 41	PRONTO-250-PLUS

**Example**

Command: *GCR	Answer: Range: 10<CR><LF>
---------------	---------------------------

05 - Set Autoscale

This command is used to force the display into autoscale.

Command	Parameters	Answer	Model Available
SAS	1: On 0: Off		PRONTO-250-PLUS

06 - Get Autoscale

This command returns whether or not the autoscale option is activated.

Command	Parameters	Answer	Model Available
GAS	None	1: On 0: Off	PRONTO-250-PLUS

**Example**

Command: *GAS	Answer: AutoScale: 1<CR><LF>
---------------	------------------------------

07 - Display Valid Scale

This command is used to display all the valid scales that the connected head supports. The scales are displayed in scale index. Please refer to the Set Scale section for the table correspondence.

Command	Parameters	Answer	Model Available
DVS	None	The valid scale index.	PRONTO-250-PLUS

The following example is for a PRONTO-250-PLUS, which can have the following scales:

1 J  
3 J  
10 J  
30 J  
100 J  
300 J

**Example**

Command: *DVS	Answer: [24]: 1.000<CR><LF> [25]: 3.000<CR><LF> [26]: 10.00<CR><LF> [27]: 30.00<CR><LF> [28]: 100.0<CR><LF> [29]: 300.0<CR><LF>
---------------	--

08 - Set Trigger Level

This command sets the internal trigger level when using the device in energy mode.

Command	Parameters	Answer	Model Available
STL	Trigger Level (in percentage) must be 4 numerical values		PRONTO-250-PLUS

**Default:** 2%

The value should be set between 0.1 and 99.9.

**Example**

Command: *STL15.4 (15.4%) *STL00.2 (0.2%)	Answer:
--	---------

09 - Get Trigger Level

This command returns the trigger level in %. The value is between 0.1% and 99.9%.  
*This is for PRONTO in Energy mode only.*

Command	Parameters	Answer	Model Available
GTL	None	Returns the trigger level in %.	PRONTO-250-PLUS

**Example**

Command: *GTL	Answer: 2.0<CR><LF>
---------------	---------------------

10 – Get Measure Mode Display

This command returns the PRONTO's measurement mode. Depending on the type of PRONTO, it can be Continuous Power Mode in W (CWP), Single Shot Power Mode in W (SSP) or Single Shot Energy Mode in J (SSE).

Command	Parameters	Answer	Model Available
GMD	None	CWP = 0 SSP = 3 SSE = 2	All

**Example**

Command: *GMD	Answer: Mode: 0<CR><LF>
---------------	-------------------------

11 – Control LCD

This command is used to control the LCD by turning it ON and OFF. The PRONTO is still working even if the LCD is off.

Command	Parameters	Answer	Model Available
LCD	1: On 0: Off		All

**Example**

Command: *LCD1	Answer:
----------------	---------

### 5.5.2. Data Acquisition

#### 12 - Query Current Value

This command is used to query the value that is currently being displayed by the monitor. The value is displayed in watts or in joules.

Command	Parameters	Answer	Model Available
CVU	None	Data in ASCII (Scientific notation)	All



#### Examples

For example, a 506.601 watts reading and a -12.25631 milliwatts reading would be displayed like this:

Command: *CVU	Answer: +5.066010e+02<CR><LF>
Command: *CVU	Answer: -1.225631e-02<CR><LF>

#### 13 - Send Continuous Transmission of Data

This command is used to send data to the serial port at a frequency of 6.8Hz in CWP mode.

Command	Parameters	Answer	Model Available
CAU	None	Data in ASCII (Scientific notation)	All



#### Examples

For example, with a PRONTO in CWP, a reading around 5.0 watts would be displayed like this until the command \*CSU is sent:

Command: *CAU	Answer : +5.066010e+00<CR><LF> +5.066012e+00<CR><LF> +5.066014e+00<CR><LF> +5.066022e+00<CR><LF> +5.066032e+00<CR><LF> +5.066042e+00<CR><LF> ...
---------------	---

14 - Stop the CAU Command

This command is used to stop the real time transfer enabled by the CAU command.

Command	Parameters	Answer	Model Available
CSU	None		All

15 - Query New Value Ready

This command is used to check whether a new value is available from the device. Though optional, its use is recommended when used with single pulse operation

Command	Parameters	Answer	Model Available
NVU	None	New Data Available or New Data Not Available	All

**Example**

Command: *NVU	Answer: New Data Not Available <CR><LF>
---------------	---

16- Send Temperature of Detector

This command is used to send temperature's data to the serial port according to the data sampling rate. To use the TMP1 command, the CAU command needs to be activated. Using TMP1 command will send temperature at the same rate as CAU command. The first number is the measured power and the second number is the measured temperature. TMP0 will send a single value of the detector temperature.

Command	Parameters	Answer	Model Available
TMP	0,1	Data in ASCII	All

**Examples**

For example, with a wattmeter, a reading around 5.0 watts would be displayed like this until the command \*CSU is sent:

Command: *CAU and *TMP1	Answer : +5.066010e+00, 23.1 <CR><LF> +5.066012e+00, 23.2 <CR><LF> +5.066014e+00, 23.2 <CR><LF>
Command: *TMP0	Answer : 23.0 <CR><LF>

### 5.5.3. Setup

#### 17 - Set Personal Wavelength Correction in nm

This command is used to specify the wavelength in nm being used on the detector. The internal memory in the detector contains measured spectral data for a wide range of wavelengths. A valid value is set between the lowest and highest wavelengths supported by the device, and it should not be a floating point value. The input parameter must have 5 digits. If the desired wavelength does not have 5 digits you must enter a zero-padded number. For example, to set the wavelength at 514 nm, you must enter 00514 or 514.0.

Specifying zero as a wavelength or providing an out-of-bound value as a parameter cancels the command.

Command	Parameters	Answer	Model Available
PWC	Wavelength		All

**Default:** Calibration wavelength, (typically 1064 nm, varies with the detector model)



#### Example

The following example sets the wavelength to 1550 nm.

Command: *PWC01550	Answer:
--------------------	---------

#### 18 - Set Personal Wavelength Correction in microns

This command is used to specify the wavelength in microns. The internal memory in the detector contains measured spectral data for a wide range of wavelengths. A valid value is set between the lowest and highest wavelengths supported by the device. The input parameter must have 5 digits and can be a floating point value. If the desired wavelength does not have 5 digits you must enter a zero-padded number. For example, to set the wavelength at 10.60 micron, you must enter 010.6.

Specifying zero as a wavelength or providing an out-of-bound value as a parameter cancels the command.

Command	Parameters	Answer	Model Available
PWM	Wavelength		All

**Default:** Calibration wavelength, (typically 1064 nm, varies with the detector model)



#### Example

The following example sets the wavelength to 2.5 micron (2500 nm).

Command: *PWM02.50	Answer:
--------------------	---------

19 - Get Wavelength

This command returns the wavelength in nm.

Command	Parameters	Answer	Model Available
GWL	None	Returns the wavelength in nm	All

**Example**

Command: *GWL	Answer: PWC: 1064<CR><LF>
---------------	---------------------------

**5.5.4. Control**20 - Set Anticipation

This command is used to enable or disable the anticipation processing when the device is reading. The anticipation is a software-based acceleration algorithm that provides faster readings using the detector's calibration.

Command	Parameters	Answer	Model Available
ANT	1: On 0: Off		All

**Default:** On

**Example**

The following example sets the anticipation On.

Command: *ANT1	Answer:
----------------	---------

21 - Get Anticipation Status

This command returns the anticipation status. If the anticipation is not available, it will always be at "off".

Command	Parameters	Answer	Model Available
GAN	None	1: On 0: Off	All

**Example**

Command: *GAN	Answer: Anticipation: 0<CR><LF>
---------------	---------------------------------



22 - Set Zero Offset

This command subtracts the current value from all future measurements the moment the command is issued to set a new zero point.

Command	Parameters	Answer	Model Available
SOU	None	Autoscale: Please Wait... Done!  Fixed scale:	PRONTO-250-PLUS

**Example**

Command: *SOU	Answer: Please Wait... <CR><LF> Done! <CR><LF>
---------------	---

23 - Clear Zero Offset

This command undoes the Zero Offset command to set the zero point at zero (cancel the SOU command).

Command	Parameters	Answer	Model Available
COU	None		PRONTO-250-PLUS

24 - Get Zero Offset

This command returns whether the zero offset has been activated or not.

Command	Parameters	Answer	Model Available
GZO	None	1: On 0: Off	PRONTO-250-PLUS

**Example**

Command: *GZO	Answer: Zero: 0<CR><LF>
---------------	-------------------------

25 - Set User Multiplier

This command is used to set the value of the multiplier between 0.5 and 2.5.

Command	Parameters	Answer	Model Available
MUL	8-character numerical value		All

**Default:** 1

**Example**

The following example sets multiplier = 2.5

Command: *MUL000002.5 Or *MUL2.500000	Answer:
---	---------

26 - Get User Multiplier

This command returns the multiplier value.

Command	Parameters	Answer	Model Available
GUM	None	Current multiplier value	All

**Example**

Command: *GUM	Answer : User Multiplier: 2.5000000E+00<CR><LF>
---------------	--

27 - Set User Offset

This command is used to set the value of the offset.

Command	Parameters	Answer	Model Available
OFF	8-character numerical value		All

**Default:** 0

**Example**

The following example sets the offset to 1.5 watts or 1.5 joules.

Command: *OFF1.500000 or *OFF1.500e+0	Answer:
---	---------

The other option available is the Zero Offset. The Zero Offset operation is done first, before the User Multipliers and Offsets

28 - Get User Offset

This command returns the offset value.

Command	Parameters	Answer	Model Available
GUO	None	Current offset value	All



**Example**

Command: *GUO	Answer : User Offset: 1.5000000E+00<CR><LF>
---------------	--

29 – Short touch on button

This command is used to simulate a short touch on any button shown on the display and the home button. Each button is numbered from 1 to 7 and they are ordered like a book, from the left to right and top to bottom. The 0 number is reserved to control the home button.



Note: This command is only supported for the following example.

Command	Parameters	Answer	Model Available
PRS	0,1,2,3,4,5,6,7	ACK or NACK	All



**Example**



The following example shows how to get the device ready for a measurement and how to save data.

Command: *PRS1 from the main menu 	Answer :ACK 
---	---



**Example**

The following example shows how to save data.

Command: *PRS2 from the main menu 	Answer : ACK 
---	--



Example

The next example shows how to change the mode using the following sequence (For PRONTO-250-PLUS).

Command:  
\*PRS0 from the main menu

Answer :ACK

\*PRS4 from the main menu

Answer : ACK

Then use PRS with one of the following number

PRS1=SSP  
PRS2=CWP  
PRS3=SSE

30 – Long touch on button

This command is used to simulate a long touch of 3 seconds on any button shown on the display and on the home button. Each button is numbered from 1 to 7 and they are ordered like a book, from the left to right and top to bottom. The 0 number is reserved to control the home button.

Note: This command is only supported for the following example

Command	Parameters	Answer	Model Available
PRL	0,1,2,3,4,5,6,7	ACK or NACK	All



Example

Command: *PRL0 from the any menu	Answer : ACK (PRONTO shut down)
*PRL2 from the main menu	ACK (Erase data, see section 2.6.3)

### 5.5.5. Instrument and Detector Information

#### 31 - Query Version

This command is used to query the device to get information about the firmware version and the device type.

Command	Parameters	Answer	Model Available
VER	None	Version and device type	All



#### Example

Command: *VER	Answer: Pronto Plus Version 1.00.04 <CR><LF>
---------------	---

#### 32 - Query Status

This command is used to query the device to get information about the following characteristics:

- Measure Mode
- Maximum, minimum and current scale
- Maximum, minimum and current wavelength with and without attenuation
- Attenuator availability and status
- Detector model
- Detector serial number

Command	Parameters	Answer	Model Available
STS	None	A hexadecimal structure described in the table below.	All

The first byte represents the validity of the structure: 0 represents a valid line while 1 is the end of the structure. The next 4 bytes represent the address line and the last 4 bytes are the actual value. The values are written on 32 bits, which means that all the values are written on two lines. The first line represents the LSB and the second line represents the MSB.

The following table shows the output WITH a XLP12-3S-H2-INT-D0 s/n 199672 (Also works with the PRONTO).

Hexadecimal Structure			Converted Value	Definition
Valid	Address	Value		
:0	0000	0003	3	Reserved
:0	0001	0000	0	Reserved
:0	0002	0003	3	Reserved
:0	0003	0000	0	Reserved
:0	0004	0000	0	Measure Mode LSB
:0	0005	0000	0	Measure Mode MSB
:0	0006	0015	21	Current scale LSB (refer to scale index *SCS)
:0	0007	0000	0	Current scale MSB (refer to scale index *SCS)
:0	0008	0019	25	Maximum scale LSB (refer to scale index *SCS)
:0	0009	0000	0	Maximum scale MSB (refer to scale index *SCS)
:0	000A	0011	17	Minimum scale LSB (refer to scale index *SCS)
:0	000B	0000	0	Minimum scale MSB (refer to scale index *SCS)
:0	000C	0428	1064	Current wavelength LSB (nm)
:0	000D	0000	0	Current wavelength MSB (nm)
:0	000E	2968	10600	Maximum wavelength LSB (nm)

:0	000F	0000	0	Maximum wavelength MSB (nm)
:0	0010	00C1	193	Minimum wavelength LSB (nm)
:0	0011	0000	0	Minimum wavelength MSB (nm)
:0	0012	0001	1	Is Attenuator available LSB (1= yes 0 = no)
:0	0013	0000	0	Is Attenuator available MSB (1= yes 0 = no)
:0	0014	0000	0	Is Attenuator on LSB (1= yes 0 = no)
:0	0015	0000	0	Is Attenuator on MSB (1= yes 0 = no)
:0	0016	2968	10600	Maximum wavelength with attenuation LSB (nm)
:0	0017	0000	0	Maximum wavelength with attenuation MSB (nm)
:0	0018	00C1	193	Minimum wavelength with attenuation LSB (nm)
:0	0019	0000	0	Minimum wavelength with attenuation MSB (nm)
:0	001A	4C 58	X L	Detector name (You must convert the hexadecimal values in ASCII characters)
:0	001B	31 50	P 1	
:0	001C	2D 32	2 -	
:0	001D	53 33	3 S	
:0	001E	- H		
:0	001F	2 -		
:0	0020	D 0		
:0	0021			0000 = Null termination character
:0	0022			The rest of the characters aren't valid until line 002A
:0	0023			
:0	0024			
:0	0025	40 03	@	
:0	0026	00 1A		
:0	0027	00 00		
:0	0028	E1 20	á	
:0	0029	00 3A	:	
:0	002A	39 31	1 9	Detector serial number (You must convert the hexadecimal values in ASCII characters)
:0	002B	36 39	9 6	
:0	002C	32 37	7 2	
:0	002D	00 00		0000 = Null termination character
:1	0000	00 00		End of structure

### 33 - Query Extended Status

This command is used to query the device to get information about the following characteristics:

Measure Mode  
Maximum, minimum and current scale  
Maximum, minimum and current wavelength with and without attenuation  
Attenuator availability and status  
Detector model  
Detector serial number  
Trigger level (0.001 to 0.999)  
Autoscale mode  
Anticipation mode  
Zero offset mode  
User multiplier  
User offset

Command	Parameters	Answer	Model Available
ST2	None	A hexadecimal structure described in the table below.	All

The first byte represents the validity of the structure: 0 represents a valid line while 1 is the end of the structure. The next 4 bytes represent the address line and the last 4 bytes are the actual value. The values are written on 32 bits, which means that all the values are written on two lines. The first line represents the LSB and the second line represents the MSB.

The following table shows the output WITH a XLP12-3S-H2-INT-D0 s/n 199672 (Also works with the PRONTO).

Hexadecimal Structure			Converted Value	Definition
Valid	Address	Value		
:0	0000	3	3	Reserved
:0	0001	0	0	Reserved
:0	0002	3	3	Reserved
:0	0003	0	0	Reserved
:0	0004	0	0	Measure Mode LSB
:0	0005	0	0	Measure Mode MSB
:0	0006	11	17	Current scale LSB (refer to scale index *SCS)
:0	7	0	0	Current scale MSB (refer to scale index *SCS)
:0	0008	19	25	Maximum scale LSB (refer to scale index *SCS)
:0	0009	0	0	Maximum scale MSB (refer to scale index *SCS)
:0	000A	11	17	Minimum scale LSB (refer to scale index *SCS)
:0	000B	0	0	Minimum scale MSB (refer to scale index *SCS)
:0	000C	428	1064	Current wavelength LSB (nm)
:0	000D	0	0	Current wavelength MSB (nm)
:0	000E	2968	10600	Maximum wavelength LSB (nm)
:0	000F	0	0	Maximum wavelength MSB (nm)
:0	0010	00C1	193	Minimum wavelength LSB (nm)
:0	0011	0	0	Minimum wavelength MSB (nm)
:0	0012	1	1	Is Attenuator available LSB (1= yes 0 = no)
:0	0013	0	0	Is Attenuator available MSB (1= yes 0 = no)
:0	0014	0	0	Is Attenuator on LSB (1= yes 0 = no)
:0	0015	0	0	Is Attenuator on MSB (1= yes 0 = no)
:0	0016	2968	10600	Maximum wavelength with attenuation LSB (nm)
:0	0017	0	0	Maximum wavelength with attenuation MSB (nm)
:0	0018	00C1	193	Minimum wavelength with attenuation LSB (nm)
:0	0019	0	0	Minimum wavelength with attenuation MSB (nm)
:0	001A	4C 58	X L	Detector name (You must convert the hexadecimal values in ASCII characters)
:0	001B	31 50	P 1	
:0	001C	2D 32	2 -	
:0	001D	53 33	3 S	
:0	001E	48 2D	- H	
:0	001F	2D 32	2 -	
:0	0020	30 44	D 0	
:0	0021	0 0		0000 = Null termination character
:0	0022	0 0		The rest of the characters aren't valid until line 002A
:0	0023	0 0		
:0	0024	1F 0		
:0	0025	40 3	@	
:0	0026	0 1A		
:0	0027	0 0		
:0	0028	E1 20	Å	
:0	0029	0 3A	:	
:0	002A	39 31	1 9	Detector name (You must convert the hexadecimal values in ASCII characters)
:0	002B	36 39	9 6	
:0	002C	32 37	7 2	
:0	002D	0 0		0000 = Null termination character
:0	002E	D70A	0.0200	Trigger Level LSB (between 0.001 and 0.999)
:0	002F	3CA3		Trigger Level MSB (between 0.001 and 0.999)
:0	0030	0001	1	Is autoscale mode on? LSB
:0	0031	0000	0	Is autoscale mode on? MSB
:0	0032	0000	0	Is anticipation on? LSB
:0	0033	0000	0	Is anticipation on? MSB
:0	0034	0000	0	Is zero offset on? LSB

:0	0035	0000	0	Is zero offset on? MSB
:0	0036	0000	1.0000	Correction Multiplier LSB
:0	0037	3F80		Correction Multiplier MSB
:0	0038	0000	0.0000	Correction Offset LSB
:0	0039	0000		Correction Offset MSB
:1	0000	0000	0	End of structure

### 34 – Return Instrument ID

This command is used to get the information about the device type.

Command	Parameters	Answer	Model Available
IDN	None	Device type	All



#### **Example**

Command: *IDN	Answer: Pronto Plus <CR><LF>
---------------	------------------------------

### 35 – Return Instrument Firmware Version

This command is used to get the firmware version of the device.

Command	Parameters	Answer	Model Available
GSV	None	Version	All



#### **Example**

Command: *GSV	Answer: 1.00.04 <CR><LF>
---------------	--------------------------

### 36 – Return Instrument Global information

This command is used to get general information about the device. This information is included in the previous commands. The information is: the identification number of the firmware, the device model and the firmware version.

Command	Parameters	Answer	Model Available
GFW	None	Version and device type	All



#### **Example**

Command: *GFW	Answer: 104233, Pronto Plus 1.00.04-RC5 <CR><LF>
---------------	---



37 – Query Battery State

This command is used to get the battery power in percentage.

Command	Parameters	Answer	Model Available
QSO	None	Number in percentage	All

**Example**

Command: *QSO	Answer: 98 <CR><LF>
---------------	---------------------



**5.5.6. Error Messages**

#	Error	Comment
1	<b>Command Error. Command not recognized.</b>	Command is invalid.
2	<b>Command Error. Command must start with '*'</b>	All text commands must begin with a trig character (*).

## 6. USB DRIVER INSTALLATION

PRONTO USB drivers will install a virtual COM port on your PC. Please download the USB driver at: <https://gentec-eo.com/downloads>.

1. Do not connect the PRONTO to your computer
2. Follow the installation steps until you have the message PRONTO ready to use.

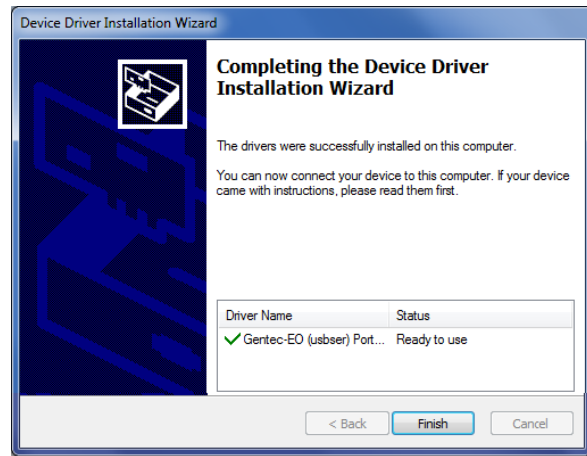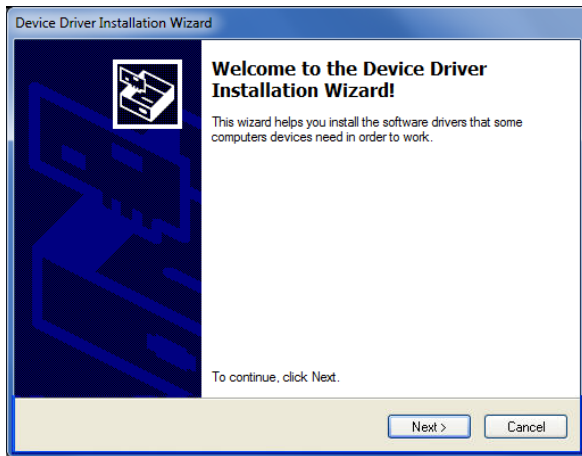

3. You can now connect the PRONTO and install the software.

## 7. MAINTENANCE

### 7.1. FREE FIRMWARE UPGRADE

As new and improved versions of the device's firmware are created, it is in your best interest to update your PRONTO. The latest device firmware can be downloaded from the Gentec-EO website at <https://gentec-eo.com/downloads>. Go to the **Downloads** section. Find the file that corresponds to your PRONTO and follow our simple, easy to use instructions.

## 8. DECLARATION OF CONFORMITY

Application of Council Directive(s): 2014/30/EU EMC Directive

Manufacturer's Name: Gentec Electro Optics, Inc.  
 Manufacturer's Address: 445 St-Jean Baptiste, suite 160  
 (Québec), Canada G2E 5N7

European Representative Name: Laser Components S.A.S.  
 Representative's Address: 45 bis Route des Gardes  
 92190 Meudon (France)

Type of Equipment: Optical Power Monitor  
 Model No.: PRONTO Series  
 Year of test & manufacture: 2015



### Standard(s) to which Conformity is declared:

#### Emissions:

Product Standard	Test Standard	Description
EN 61326-1_Ed2:2013 (IEC 61326-1_Ed2:2012)	CISPR 11:+A1:2010 Class A	Radiated Emissions
EN 61326-1_Ed2:2013 (IEC 61326-1_Ed2:2012)	Radiated Emissions FCC part 15 2013) subpart B	Radiated Emissions

#### Immunity:

Product Standard	Test Standard	Description	Performance Criteria
EN 61326-1_Ed2:2013 (IEC 61326-1_Ed2:2012)	IEC 61000-4-2:2008Ed.2	Electrostatic Discharge Immunity	Criteria B
EN 61326-1_Ed2:2013 (IEC 61326-1_Ed2:2012)	IEC 61000-4-3:2006+ A1:2007+A2:2010	RF Conducted Immunity	Criteria A

I, the undersigned, hereby declare that the equipment specified above conforms to the above Directive(s) and Standard(s).

Place: Québec (Québec)

Date : 7 July, 2015

## 9. UKCA DECLARATION OF CONFORMITY

Application of Council Directive(s): 2014/30/EU EMC Directive

Manufacturer's Name: Gentec Electro Optics, Inc.  
 Manufacturer's Address: 445 St-Jean Baptiste, suite 160  
 (Québec), Canada G2E 5N7

European Representative Name: Laser Components S.A.S.  
 Representative's Address: 45 bis Route des Gardes  
 92190 Meudon (France)

Type of Equipment: Optical Power Monitor  
 Model No.: PRONTO Series  
 Year of test & manufacture: 2015



### Standard(s) to which Conformity is declared:

#### Emissions:

Product Standard	Test Standard	Description
EN 61326-1_Ed2:2013 (IEC 61326-1_Ed2:2012)	CISPR 11:+A1:2010 Class A	Radiated Emissions
EN 61326-1_Ed2:2013 (IEC 61326-1_Ed2:2012)	Radiated Emissions FCC part 15 2013) subpart B	Radiated Emissions

#### Immunity:

Product Standard	Test Standard	Description	Performance Criteria
EN 61326-1_Ed2:2013 (IEC 61326-1_Ed2:2012)	IEC 61000-4-2:2008Ed.2	Electrostatic Discharge Immunity	Criteria B
EN 61326-1_Ed2:2013 (IEC 61326-1_Ed2:2012)	IEC 61000-4-3:2006+ A1:2007+A2:2010	RF Conducted Immunity	Criteria A

I, the undersigned, hereby declare that the equipment specified above conforms to the above Directive(s) and Standard(s).

Place: Québec (Québec)

Date : December 01, 2021

## APPENDIX A: WEEE DIRECTIVE

### Recycling and separation procedure for WEEE directive 2012/19/EU

This section is used by the recycling center when the detector reaches the end of its life. Removing the insulation or troubling the inside of the monitor will void the detector warranty.

The complete Detector contains:

- 1 Detector
- 1 Cable
- 1 Calibration certificate

### Separation

Paper: Certificate

Printed Circuit Board: Inside the Detector

Aluminum: Detector Casing

Plastic: Parts Inside the Detector


Li-ion cell: Battery

## APPENDIX B: INSTALLING THE PRONTODATATRANSFER SOFTWARE


1. Download and install the driver from our website.
2. Download the software file from our website.
3. Double-click the EXE file to start the installer.




4. Choose the folder for the installation and click Next until the installation is complete. Then click Close.

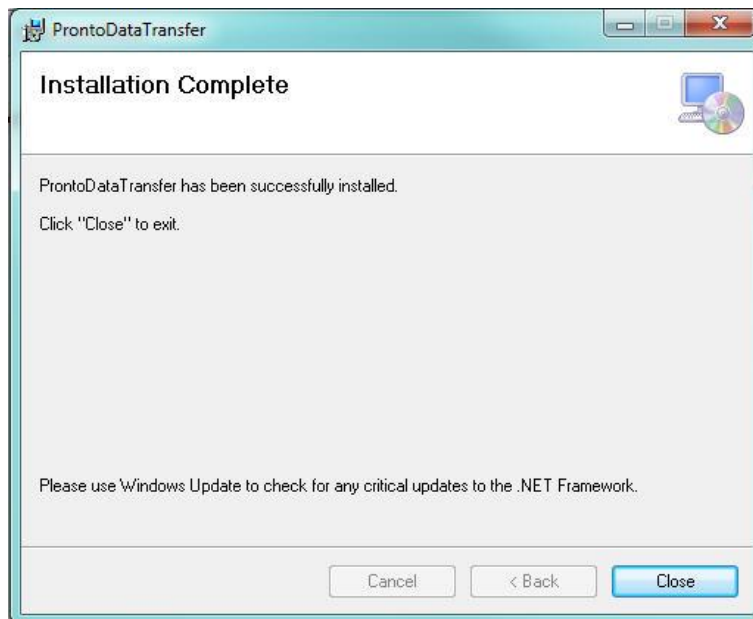


5. Once the software is installed, locate it on your computer and start the program by clicking the

PRONTODataTransfer icon .


6. A download window appears with a **Download File** button.

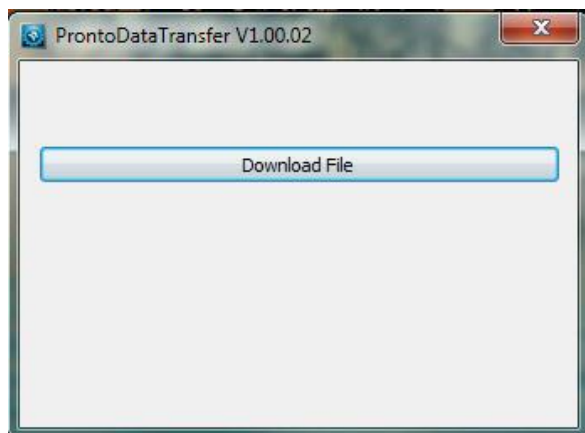

7. Click the **Download File** button and select a known folder on your computer where you will save the data. Don't forget to also enter a file name. The file format is TXT.

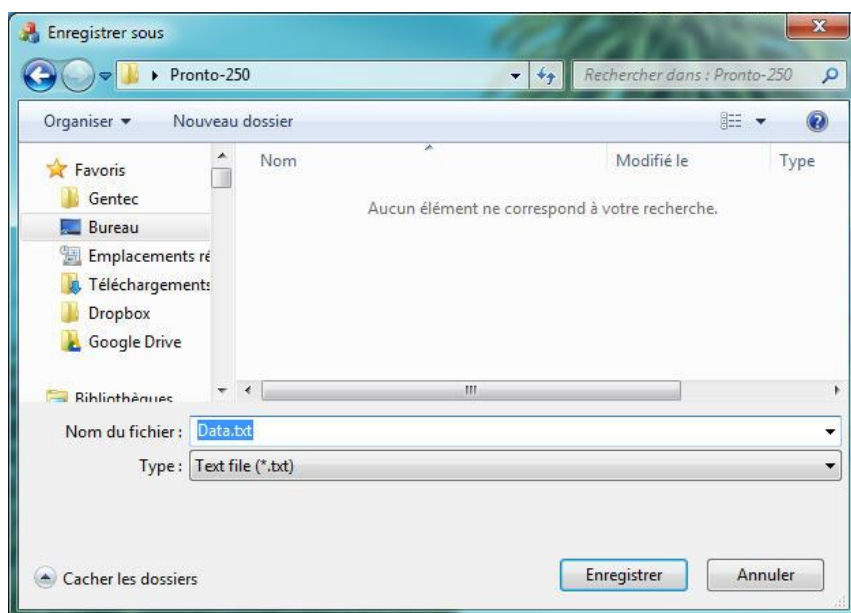

8. You can now open and analyze the data in your preferred data analysis software.

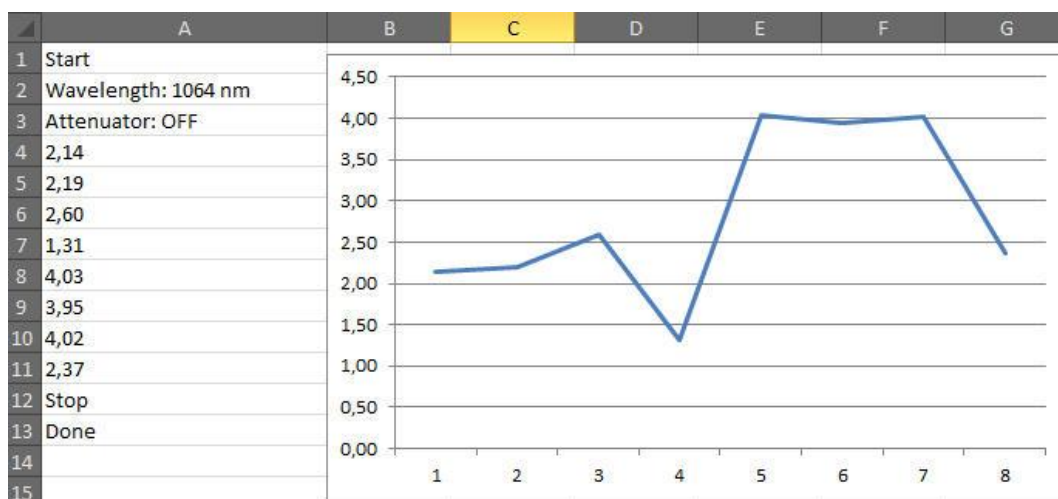

## APPENDIX C: UPDATING THE PRONTO FIRMWARE


1. Before updating the firmware, transfer any saved data using the PRONTODataTransfer software.
2. Download the executable from our website.
3. Plug your PRONTO to your computer using a USB cable.




### Warning

Do not unplug the PRONTO or turn of your computer while you are updating your device.

4. Double-click the EXE file to start the updater. Click Yes to begin the update.

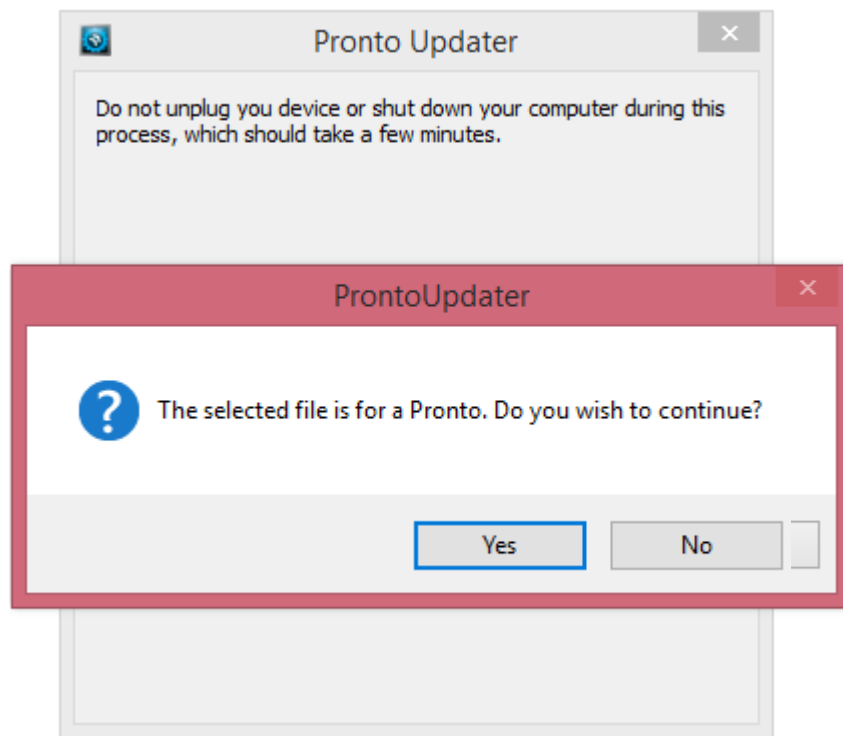

5. Once the update is done, you can unplug your device and continue using your PRONTO as usual.

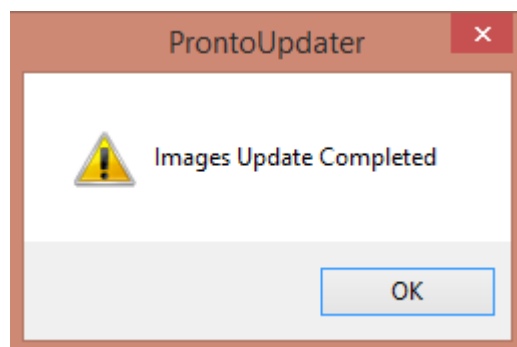





**gentec-eo**  
PARTNERS for ACCURACY

[WWW.GENTEC-EO.COM](http://WWW.GENTEC-EO.COM)

## LEADER IN LASER BEAM MEASUREMENT SINCE 1972



POWER & ENERGY METERS



BEAM PROFILING



THZ MEASUREMENT

### CANADA

445 St-Jean-Baptiste, Suite 160  
Quebec, QC, G2E 5N7  
CANADA

T (418) 651-8003  
F (418) 651-1174

[info@gentec-eo.com](mailto:info@gentec-eo.com)

### UNITED STATES

5825 Jean Road Center  
Lake Oswego, OR, 97035  
USA

T (503) 697-1870  
F (503) 697-0633

[info@gentec-eo.com](mailto:info@gentec-eo.com)

### JAPAN

Office No. 101, EXL111 building,  
Takinogawa, Kita-ku, Tokyo  
114-0023, JAPAN

T +81-3-5972-1290  
F +81-3-5972-1291

[info@gentec-eo.com](mailto:info@gentec-eo.com)

### CALIBRATION CENTERS

- 445 St-Jean-Baptiste, Suite 160  
Quebec, QC, G2E 5N7, CANADA
- Werner von Siemens Str. 15  
82140 Olching, GERMANY
- Office No. 101, EXL111 building,  
Takinogawa, Kita-ku, Tokyo  
114-0023, JAPAN